

SAMPLING INSTRUCTIONS

The following sampling instructions are used for collecting water or groundwater samples for plate count and/or most probable number (MPN) analysis. The recommended sampling container is two unpreserved 40mL VOA vials for each sampling location. Microbial Insights, Inc. can provide the proper sampling supplies upon request.

Once the proper sampling vial is obtained be sure not to contaminate the inside of the sample vial with skin, dirt or any form of debris (this helps to ensure the accuracy of the data results). Wearing latex gloves (or similar) is recommended when sampling.

Please fill each of the 40mL VOA vials with the sampling water leaving no headspace in the vial.

* Note: It is important to collect as close to the required amounts as possible to ensure the ability to properly conduct the analysis requested. Hold time for this analysis is 24 – 48 hours.

To Submit Sample:

1. Once the required amount of groundwater has been collected into the proper sampling container, seal the container tightly with a screw cap lid.
2. Properly affix a label with the sample name, date and analysis.
3. Be sure to fill out the Chain of Custody (COC) form correctly and accurately and ship it along with the samples. A COC form is required for QA/QC purposes.
4. Once the vials have been correctly labeled, place them in the designated cooler. Be sure to fill the remaining space in the cooler with blue ice or regular ice that has been double bagged in Ziploc bags to keep the entire shipment around 4°C, especially during the summer months.
5. All paperwork to be sent with the samples should be placed within a waterproof pouch or Ziploc bag and placed on top of the samples or affixed to the inside lid of the cooler.
6. Seal the cooler lid with a strong packaging tape.

SHIPPING INSTRUCTIONS

Packaging Samples:

1. Samples should be shipped in a cooler with ice or blue ice for next day delivery. If regular ice is used, the ice should be double bagged.
2. A chain of custody form must be included with each shipment of samples. Access our chain of custody at www.microbe.com.

Shipment for Weekday Delivery:

Samples for weekday delivery should be shipped to:

Sample Custodian
Microbial Insights, Inc.
10515 Research Drive
Knoxville, TN 37932
(865) 573-8188

Shipment for Saturday Delivery:

Note: Microbial Insights, Inc is **closed** on Sunday, however we can receive samples on Saturday. Please contact us prior to shipping if the delivery of the samples is going to be on a Saturday.

Samples for **Saturday delivery** should be shipped to:

Microbial Insights, Inc.
FedEx Drop Location
10601 Murdock Drive
Knoxville, TN 37932
(865) 300-8053